


Instructions & Sample Pit Survey Diagram Showing Locations

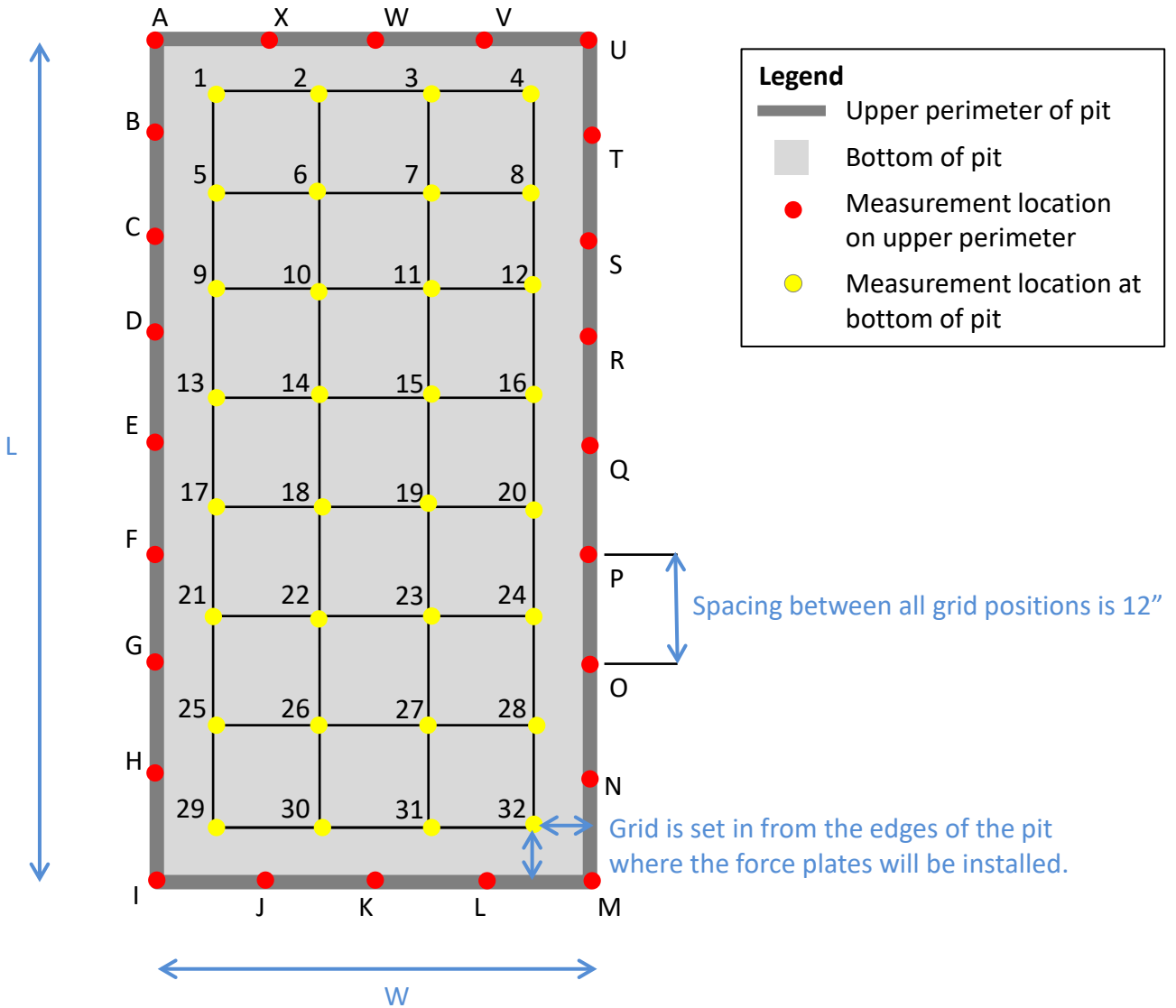

Position a laser level on the surrounding floor. Measure the depth from the plane to the floor along the upper perimeter of the pit on 12" increments and at every corner. Identify the high spot.

Without moving the laser level measure the depth to the bottom of the pit on a 12" grid. Focus on the area in which the plates will be located, as shown

Submit all raw data to AMTI marked on a sketch or using a labeled sketch and table, as shown. This can be hand sketched and hand written

Please provide a sketch that includes the overall dimensions of your pit and the location of the grid within the pit

Sample Table for Reporting Survey Values

Enter the distance from the reference plane to the pit perimeter or base at each location

Position on Perimeter	Distance from laser level (indicate units)
A	
B	
C	
D	
E	
F	
G	
H	
I	
J	
K	
L	
M	
N	
O	
P	
Q	
R	
S	
T	
U	
V	
W	
X	

Position at base of pit	Distance from laser level (indicate units)
1	
2	
3	
4	
5	
6	
7	
8	
9	
10	
11	
12	
13	
14	
15	
16	
17	
18	
19	
20	
21	
22	
23	
24	
25	
26	
27	
28	
29	
30	
31	
32	